

GGNB Method Course

PCR:
self-made enzymes,
helpful additives
and insights into the reactions

PRACTICAL PART

16.10.2012

Koray Kirli and Steffen Frey
Cellular Logistics
Prof. Dirk Görlich
MPI for Biophysical Chemistry
Am Fassberg 11
37077 Göttinge

I. Expression

	PfuS-Polymerase	Pab-PPase	Pab-dUTPase
o/n culture	pSF421 in BL21 DE3 Star pRosetta, 0,5 l 2YT/ 300µg/ml Amp/ 34µg/ml CAM, 37°C	pSF318 in BL21 DE3 Star pRosetta, 0,5 l 2YT/ 300µg/ml Amp/ 34µg/ml CAM, 37°C	pSF323 in BL21 DE3 Star pRosetta, 0,5 l 2YT/ 2% glycerol/ 300µg/ml Amp/ 34µg/ml CAM, 37°C
Dilution	To 3,5l with prewarmed (37°C) TB/Amp/Cam final OD ₆₀₀ : 0.5-0.8 distribute in 4x 850ml in 5l flasks	To 3,5l with prewarmed (37°C) TB/Amp/Cam final OD ₆₀₀ : 0.5-0.8 distribute in 4x 850ml in 5l flasks	To 3,2l with prewarmed (37°C) 2YT/Amp/Cam + 400mM NaCl final OD ₆₀₀ : 0.3 distribute in 4x 800ml in 5l flasks
Shaking	37°C, 100rpm	37°C, 100rpm	37°C, 100rpm
Induction	at OD ₆₀₀ =3.0 add 0.2mM IPTG final	at OD ₆₀₀ =3.0 add 0.2mM IPTG final	at OD ₆₀₀ =0.8 add 1mM IPTG final
Harvest	3-4h after induction add 1mM PMSF (stock 100mM in EtOH) centrifuge 7min, 7000rpm	3-4h after induction add 1mM PMSF (stock 100mM in EtOH) centrifuge 7min, 7000rpm	5h after induction add 1mM PMSF (stock 100mM in EtOH) centrifuge 7min, 7000rpm
Resuspension	200 OD ₆₀₀ /ml in HS-buffer + 1mg/ml Lysozyme + 5mM DTT + 5mM Imidazole	200 OD ₆₀₀ /ml in LS-buffer + 1mg/ml Lysozyme + 5mM DTT + 5mM Imidazole	200 OD ₆₀₀ /ml in HS-buffer + 1mg/ml Lysozyme + 5mM DTT + 5mM Imidazole + 0.1% TX-100
Lysis	freeze in liquid N2 thaw in warm water sonification 2-3x 1min on ice	freeze in liquid N2 thaw in warm water sonification 2-3x 1min on ice	freeze in liquid N2 thaw in warm water sonification 2-3x 1min on ice
Centrifugation	40000rpm, 60min	40000rpm, 60min	40000rpm, 60min
Storage	add 1/9 volume 2.4M sucrose freeze in aliquots of ≈10ml	add 1/9 volume 2.4M sucrose freeze in aliquots of ≈10ml	add 1/9 volume 2.4M sucrose freeze in aliquots of ≈10ml

II. Purification

	PfuS-Polymerase	Pab-PPase	Pab-dUTPase
thawing	RT water	see polymerases	see polymerases
sample "native"	5µl + 95µl SDS-buffer	see polymerases	see polymerases
heat denaturation	in 5x 2ml tubes 15min, 90°C in Thermomixer with 2ml-rack	see polymerases	see polymerases
centrifugation	5 min, 13000rpm, RT pool supernatant	see polymerases	see polymerases
sample "heated"	5µl + 95µl SDS-buffer	see polymerases	see polymerases
nickel-binding	add 1g Nickel-Silica 24% 500Å pre-equilibrated with HS-buffer, rotate on wheel 30min at RT	add 1g Nickel-Silica 24% 500Å pre-equilibrated with LS-buffer rotate on wheel 30min at RT	add 1g Nickel-Silica 24% 500Å pre-equilibrated with HS-buffer rotate on wheel 30min at RT
sample "flow through"	sediment beads take sample 5µl + 95µl SDS-buffer	see polymerases	see polymerases
wash 1	remove supernatant wash 2x with 10ml HS-buffer containing 5mM DTT and 5mM Imidazole	remove supernatant wash 2x with 10ml LS-buffer containing 5mM DTT and 5mM Imidazole	remove supernatant wash 2x with 10ml HS-buffer containing 5mM DTT and 5mM Imidazole
wash 2	wash 2x with 10ml LS-buffer containing 5mM DTT	wash 2x with 10ml LS-buffer containing 5mM DTT	wash 2x with 10ml HS-buffer containing 5mM DTT
transfer to column	transfer to plastic column	see polymerases	see polymerases
wash 3	wash 1x with 2ml LS-buffer containing 5mM DTT	see polymerases	wash 2 HS-buffer containing 5mM DTT
elute	elute with 10x 300µl LS-buffer + 5mM DTT + 400mM imidazole collect fractions between each fraction wait ≈ 2 min check protein content by amido black staining (0.5µl each)	see polymerases	elute with 10x 300µl HS-buffer + 5mM DTT + 600mM imidazole collect fractions between each fraction wait ≈ 2 min check protein content by amido black staining (0.5µl each)

	PfuS-Polymerase	Pab-PPase	Pab-dUTPase
pool	pool 5-7 fractions with highest protein content	see polymerases	see polymerases
sample "eluate"	5µl + 20µl SDS-buffer	see polymerases	see polymerases
protein determination	measure OD ₂₈₀ with nano-drop against elution buffer PfuS: 1 OD ₂₈₀ = 0.73mg/ml	measure OD ₂₈₀ with nano-drop against elution buffer 1 OD ₂₈₀ = 0.70mg/ml	measure OD ₂₈₀ with nano-drop against elution buffer 1 OD ₂₈₀ = 0.75mg/ml
dilution	dilute to 1 mg/ml with 20mM Tris-HCl pH 7.5 100mM KCl 1 mM DTT 0.1 mM EDTA 0.5 % NP40 0.5 % Tween20 100µg/ml BSA 50% Glycerol	see polymerases	see polymerases

in groups of 3-4 students: prepare mixtures in dilution buffer

	PfuS	PfuS Tripple
Volume	1ml	a lot!
Polymerase	100ng/µl PfuS	100ng/µl PfuS
Pab PPase	-	15ng/µl
Pab dUTPase	-	5ng/µl

III. Activity Assay

Questions:

Effects of helper enzymes?

Speed of Polymerase mixtures?

Test-PCRs:

Template: mixture various plasmids with different insert sizes

pSF1350/787/1387/1568/961

Primers: pQE-F /pQE-R

PCR product lengths: 0.5/1kb/1.7kb/2.4kb/3kb

PCR master-mix:

50µl 5x Phusion HF buffer

25µl dNTPs 2.5mM each

2.5µl pQE-F primer

2.5µl pQE-R primer

2.5µl template-mixture (20ng/µl of each plasmid)

168µl water

PCR reactions:

aliquot 4x 50µl (on ice)

add 0.5µl polymerase mix per reaction

1: no polymerase

2: PfuS tripple from stock

3: newly-made PfuS

4: newly-made PfuS tripple

PCR program:

1	pre-heating of block	pause (∞)	98.5°C
2	initial denaturation	1 min	98.5°C
3	denaturation	20 sec	98.5°C
4	annealing	20 sec	62°C
5	elongation	30 sec	72°C
	repeat steps 3 to 5	25 times	
6	final elongation	30 sec	72°C
7	cooling	∞	10°C

IV. Quality control:

SDS gel

boil samples 5 min

load 5µl of each sample

- | | |
|---|--------------|
| 1 | empty |
| 2 | native |
| 3 | heated |
| 4 | flow through |
| 5 | marker |
| 6 | eluate |
| 7 | empty |

run 65min at 50mA, 400V

stain with Coomassie

Agarose gel

pour 1% agarose gel containing Ethidium Bromide

to each PCR reaction add 50µl DNA loading buffer

load 10µl per slot

run for 20-30min at 180V

V. Buffer compositions:

HS-buffer:

2 M NaCl

50 mM Tris-HCl (pH 8.0)

10 mM MgCl₂

LS-buffer:

44 mM Tris-HCl (pH 7.5)

290 mM NaCl

4.4 mM MgCl₂

polymerase dilution buffer:

20mM Tris-HCl pH 7.5

100mM KCl

0.1 mM EDTA

0.5 % NP40

0.5 % Tween20

100µg/ml BSA

50% Glycerol

1 mM DTT

prepare all stocks without DTT, always add DTT freshly from 1M stock.

SDS Sample buffer (1 to 2-fold)

for 250ml

62,5 ml 0.5M Tris-HCl pH 6.8

37.5 ml 20% SDS

1.93 g DTT (solid)

85.5 g sucrose (solid)

a bit of Bromphenol-Blue

Amido black solution:

0.2% Naphtol Blue Black

2% acetic acid

Anleitung für Expression,
Zusammensetzung etc ist im Anhang
zu finden. Wir verwenden den PCR-
Puffer für Phusion (HF) von NEB.
Wichtig ist, bei der PCR die
Denaturierung bei 98,5°C
durchzuführen. Die Polymerase
macht dann bis zu 6kb pro Minute.
Ich rechne routinemäßig mit 2,5-
3kb/min.